

## 12.1 Installing software on Siemens Logo!

### Preparation

- Transfer the program you want to the memory card.
- Then remove the tank sensor at the back of the EASYRO.  
System – **with level-controlled system**
- Pure water outlet of the EASYRO: Please insert a stopcock and leave it open.
- Open the water infeed.  
Note: The water infeed should always pass through a pre-filter!
- Now open the cover of the EASYRO system.
- Connect the system to the power supply.
- System displays “Check sensor” – **with level-controlled system**

### Transferring software to the control unit

First of all, the memory card must be inserted into the control unit, please proceed as follows:

- Open the card slot on the Siemens LOGO! and insert the memory card into the card slot.  
See photo



- Please close the card slot.
- Now the program must be transferred to the Siemens LOGO!; please proceed as follows:
- Press the ▼ arrow key until the date/time appears.



- Press **ESC**.
- Use the arrow keys to select “Stop” and confirm with **OK**, then select “YES” and confirm with **OK**.
- Then select the “Card” menu item and confirm with **OK**.



- Select the “Load Prog←Card” menu item and confirm with OK.



- Confirm “YES” with **OK**.



- Program is now saved from the memory card to the Siemens LOGO!
- When the process is complete, the main menu reappears on the display.
- Now switch off the power to the system and pull out the mains plug.
- Plug the mains plug back in.
- The main menu appears on the LOGO!

- Use the arrow keys to select “Start” and press , then select “YES” and confirm with .
- Close the cover.
- Changeover complete.
- Now set the correct tank height.